



LABORATORY ANALYTICAL INSTRUMENTS

pH - mV- Orp - ISE - Conducibility - TDS - Salinity - Resistivity - Temperature



MAIN FEATURES

- Simultaneous display up to 6 parameters of user's choice;
- Selectable resolution: 0.1/0.01/0.001 pH; automatic scaling for COND;
- Automatic and manual pH calibration up to 5 points;
- Automatic ISE calibration up to 5 points;
- Automatic or manual **ORP** calibration on **one point**;
- Automatic COND calibration up to 4 points and manual 1 point;
- Ultra-pure water compensation for low conductivity measurements;
- Conductivity cell with 2 and 4 rings;
- Selectable cell constant 0.1 / 1.0 / 10;
- Data recall of last calibration;
- Graphical display of **electrode condition**;
- Min-Max alarm thresholds setting for pH and ORP parameters;
- Analog view;
- Calibration due alarm;
- Data logger up to 8,000 GLP records, downloadable on PC or printer;
- Free **"DataLink+" software** for data management to PC;
- Multilanguage: Ita Eng Deu Fra Esp Por;
- USB connection for PC and RS232 for printer;



New setup menu







ADVANCED USER MANAGEMENT

The new advanced user management system includes **4 levels of access**, for 15 total profiles, with different privileges:

ADMIN - RESPONSIBLE - USER - GUEST


Through the **Datalink+ software**, available free of charge at www.xsinstruments.com, it can be viewed the **complete audit trail** with calibration information, parameter settings and active users.

GRAPHIC DISPLAY

The $8/80\ \text{PRO}$ Series features a large, backlit, $4.3\ \text{inch}\ \text{color}\ \text{display}\ \text{that}\ \text{allows}\ \text{all}\ \text{information}\ \text{to}\ \text{be}\ \text{read}\ \text{clearly}.$

Calibration points, electrode status, user and sample data, working and connection mode, logger type, date and time are displayed.

The smart interface and multilingual menu (6 selectable languages) make these tools extremely easy to use.

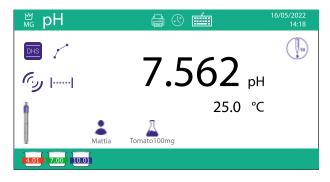

Measurement screen



EDM Ltd. โทรศัพท์ : 02 966 1590 มือถือ : 086 316 7625 Email : contact@edm.co.th LINE ID : EDM9824



The 80 PRO Series allows the use of the 4-ring conductivity cell, enabling accurate measurements over the entire scale using a single sensor.



[™] _{MG} MultiView	e 🕓 💼	16/05/2022 14:18
рН	7,000 pH	25,0 °C
mV	0,0 mV	25,0 °C
ISE	0.00 mV	25,0 °C
Cond	12.880 µS/cm	25,0 °C
TDS	0.00 mg/l	25,0 °C
SAL	0.00 ppm	25,0 °C

SETUP	É	16/05/2022 14:18
Sample ID User ID Company name Lot production Method PM name Laboratory Enviromental condition	Tomato 100mo Mattia g. Giorgio Borma 223355 305 Mattia Clinic 60,0 °C - 50%	C
to move 🔺 💎 t	o confirm 📢 👘 🕇	o exit 💮

DATA LOGGER

With four different modes of data saving, the instrument will be autonomous in monitoring the progress of analysis, while the user can focus on other operations in the laboratory:

- Manual Log;
- Automatic Log;
- Log when a Delta is reached;
- Log with Min-Max thresholds;



MAIN FEATURES

- Selectable resolution: 0.1/0.01/0.001 pH; automatic scale for COND;
- Automatic pH calibration up to 5 points and manual 2 points;
- Automatic ORP calibration on one point (475 mV);
- Automatic COND calibration up to 4 points and manual 1 point;
- Selectable cell constant 0.1 / 1 / 10;
- Data recall of last calibration;
- Graphic display of electrode condition;
- Calibration expiration alarm;
- Data logger up to 1000 GLP data with possibility of sending them to PC or printer;
- Free downloadable "DataLink+" software for data management to PC;
- Multilingual setup menu;
- USB connection for PC connection and RS232 for printer;
- Min-Max alarm thresholds setting for pH and ORP parameters;
- Numeric password for protection of calibration and saved data.
 - EDM Ltd.

MULTIVIEW

Thanks to the new **Multiview function**, up to a maximum of 6 parameters can be selected in the PC 80 PRO multiparameter to be displayed simultaneously.

The display shows in real time each parameter, measurement value and temperature.


ANALYSIS MENU

Using the Analysis menu, information regarding the sample, batch, user, supervisor, company, method, environmental conditions, etc. can be entered.

CUSTOMIZABLE TEMPLATES

Up to 4 customizable templates to tailor the measurement experience to your liking.

The templates allow you to **select which Analysis menu items to print or export to the DataLink+ software**, so you can customize your analysis reports.

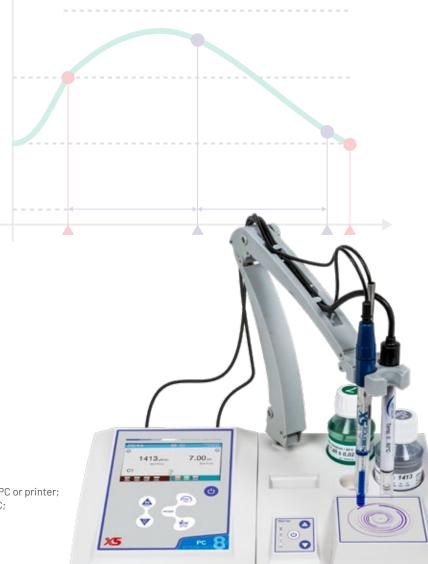

โทรศัพท์ : 02 966 1590 มือถือ : 086 316 7625 Email : contact@edm.co.th LINE ID : EDM9824

DATA SHEET	80 Pro Series	8 Pro Series	
H measuring range	-2 20	-2 16	
Resolution / accuracy	0,1 / 0,01 / 0,001 ± 0,002		
Calibration points	15		
Recognized buffers	USA - NIST - DIN - METTLER TOLEDO - MERCK - DIN/NIST + 5 CUSTOM	USA - NIST - DIN + 2 CUSTOM	
ndication of calibration points	Yes with	icons	
LP calibration	Yes with date and time		
lectrode status indication	Yes with icons and graph		
alibration expiration with alarm	Yes		
leasurement stability criteria	Medium - High - Tit (titration)		
larm settings	Min - Max		
nV measuring range (ORP) / Resolution	± 2000 0,1 / 1		
SE measuring range	0,001 19999 ppm	-	
leasuring units	Mol/L mg/L g/L	-	
alibration points	25	-	
OND measuring range	0,00 1000 mS	0,00 500 mS	
esolution	Automatic scale		
ccuracy	± 1% full scale		
OND Calibration points	14		
lecognized buffers	84 , 147 , 1413 μS , 12,88 , 111,8 mS - 1 custom value		
ndication of calibration points	Yes with icons		
emperature coefficient	Ultra-pure water - 0,00 10,00 %/°C	0,00 10,00 %/°C	
eference temperature	15 30	D°C	
LP calibration	Yes with date	and time	
alibration expiration with alarm	Yes		
ell type	4 rings - 2 rings	2 rings	
ell constant	0,475 - 0.1 - 1 - 10	0.1 - 1 - 10	
DS measuring range	0,1 mg/L 500,0 g/L	0,1 mg/L 200,0 g/L	
DS Factor/ Accuracy	0.40 1.00 / ± 1% full scale		
alinity measuring range / accuracy	0,01 100 ppt / ± 1% full scale		
Resistivity measuring range / accuracy	1 10 MΩ / ± 1% full scale	-	
emperature °C measuring range / accuracy	-30,0 130,0 °C / ±0,2 °C	-20,0 120,0 °C / ±0,2 °C	
emperature compensation	Auto e manual 0 100 °C (ΝΤC 30 ΚΩ)		
emperature calibration	Yes		
dvanced user management	4 levels / 15 profiles	-	
udit trail	Complete	-	
nalysis menu	Yes, up to 4 customizable templates	-	
lultiview	Up to 6 user-selectable parameters	pH - Cond	
assword	Alphanumeric	Numeric	
lemory	8,000 data with date and time (Man - Auto - MinMax - Delta)	1,000 data with date and time (Man - Auto)	
lisplay	Colored T	-T 4,3 "	
nputs	BNC - RCA - Multipin - Ref (banana 4mm)	BNC - RCA	
onnections	2x USB (PC - Keyboard	2x USB (PC - Keyboard) - RS232 (Printer)	
electable languages	Ita - Eng - Deu - F	ra - Esp - Por	
lagnetic stirrer	Yes, independent control		
tirring speed / Capacity	0 3000 rpm / 2 L		
^o rating	IP 54		
ower supply	AC/DC adapter 12 V / 1000 mA		
imensions (LxPxA) - Weight	165 x 195 x 68 mm - 1,05 Kg		

GIORGIO·B@RMAC





Distributed by:

SERIE_880_PR0_V1.0_05_22_ENG

EDM Ltd. EDM Ltd. [Fect c102795] YEARS โทรศัพท์ : 02 966 1590 มือถือ : 086 316 7625 Email : contact@edm.co.th LINE ID : EDM9824